



Dispenser arm kit for Frac-950

Instructions


Removing the standard mode components

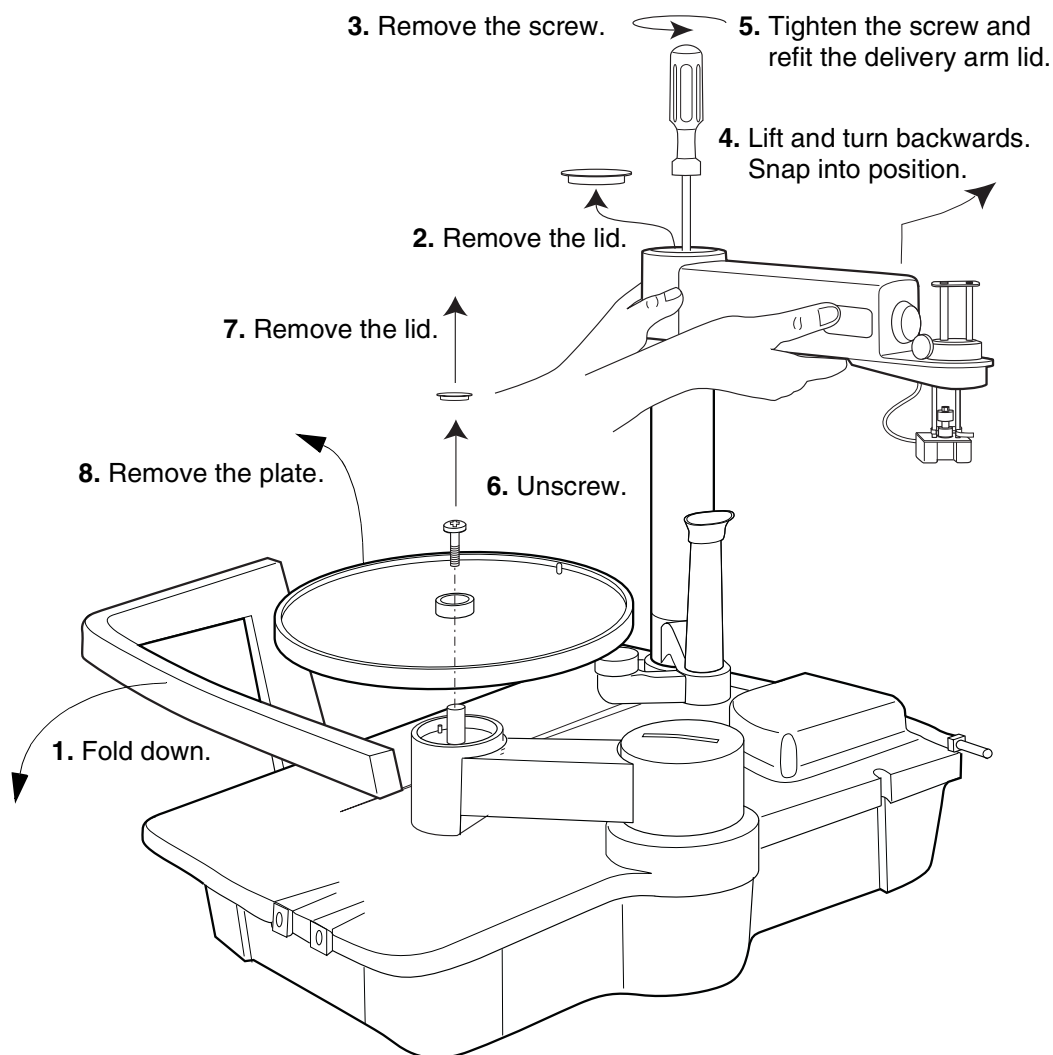

Installing the rotating shaft

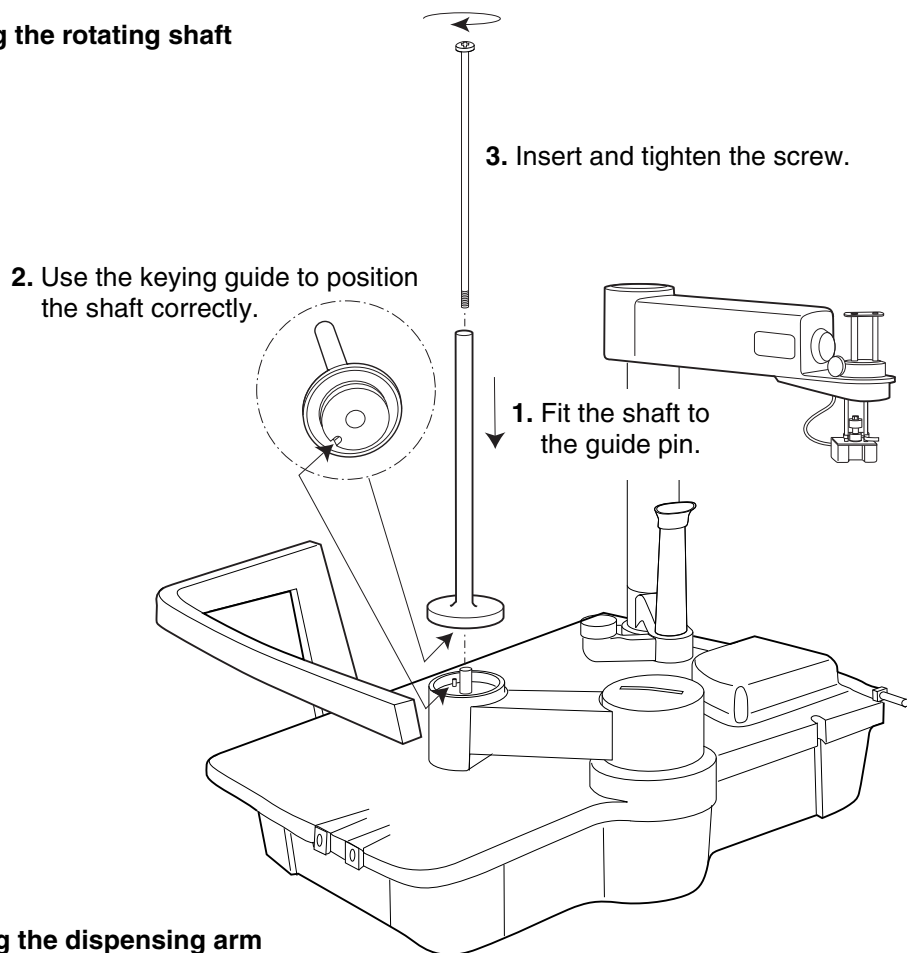

Mounting the dispensing arm

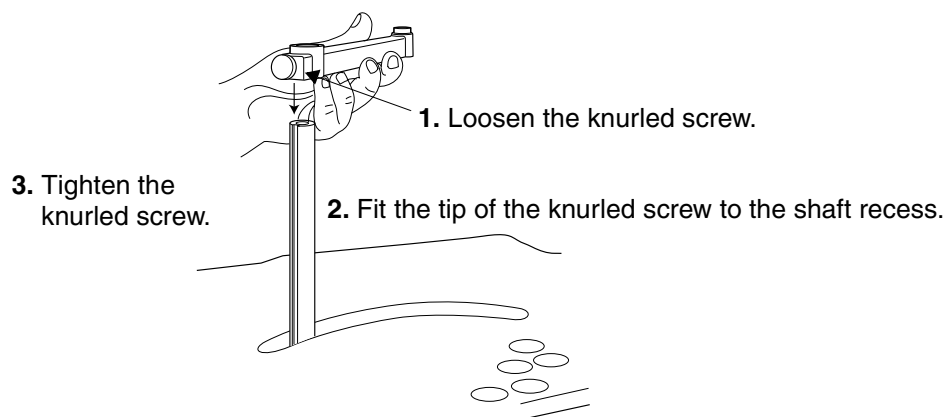


Installing the capillary

3. Thread the free end of the capillary through the fingertight connector and the swivel. Allow the capillary end to stick out 1-2 mm. Hold the swivel, with the 8 mm key supplied, and tighten the connector.

2. Fit a fingertight connector to the swivel.

4. Adjust the height of the dispenser arm to suit the height of the vessels (a distance between the arm and the vessels of 0.5-1 cm is appropriate). Tighten the knurled screw.

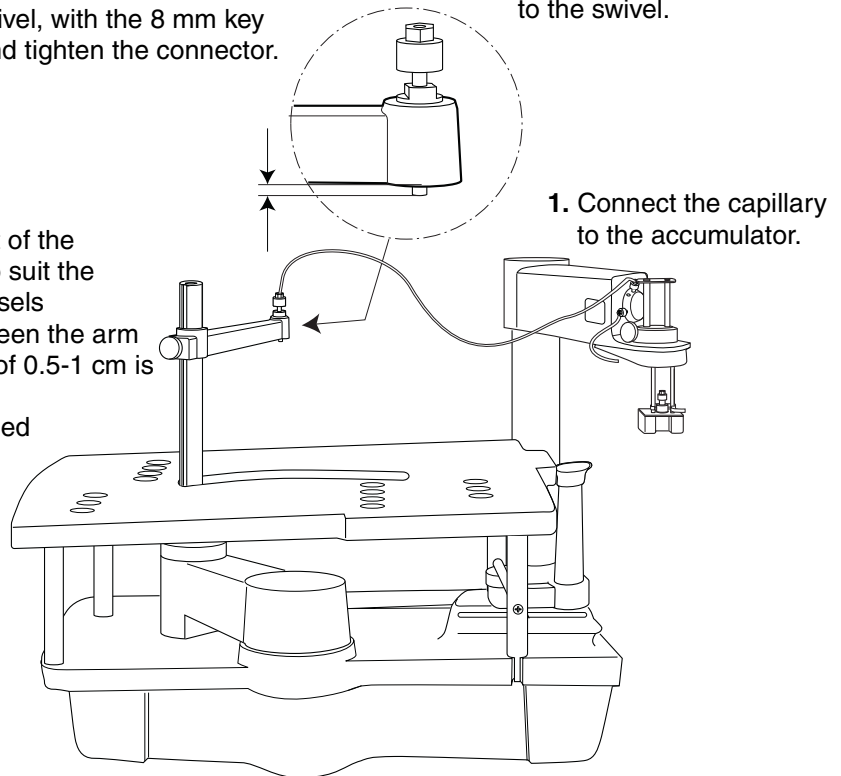

Important user information

Declaration of conformity

The **CE** symbol and corresponding declaration of conformity is valid for the instrument when it is:

- connected to other CE-marked Amersham Pharmacia Biotech instruments, or
- connected to other products recommended or described in this manual, and
- used in the same state as it was delivered from Amersham Pharmacia Biotech except for alterations described in this manual.

Terms and Conditions of Sale

All goods and services are sold subject to the terms and conditions of sale of the company within the Amersham Pharmacia Biotech group which supplies them. A copy of these terms and conditions is available on request.

Should you have any comments on this product, we will be pleased to receive them at:

Amersham Pharmacia Biotech AB

SE-751 84 Uppsala
Sweden

Trademarks

Amersham is a trademark of Nycomed Amersham plc. Pharmacia and Drop Design are trademarks of Pharmacia & Upjohn Inc.

Office Addresses

Amersham Pharmacia Biotech AB

SE-751 84 Uppsala
Sweden

Amersham Pharmacia Biotech UK Limited

Amersham Place
Little Chalfont
Buckinghamshire
England HP7 9NA

Amersham Pharmacia Biotech Inc.

800 Centennial Avenue
P.O. Box 1327
Piscataway, N.J. 08855
USA

Amersham Pharmacia Biotech Europe GmbH

Postfach 5480
D-79021 Freiburg
Germany

Amersham Pharmacia Biotech KK

Sanken Building
3-25-1 Hyakunincho, Shinjuku-ku
Tokyo 169-0073
Japan

© Copyright Amersham Pharmacia Biotech AB 2001

- All rights reserved



amersham pharmacia biotech