
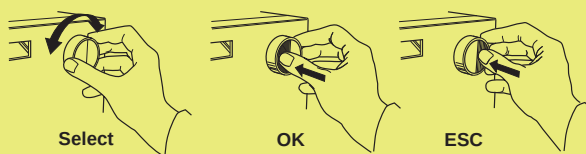


Short instructions Monitor UV-900

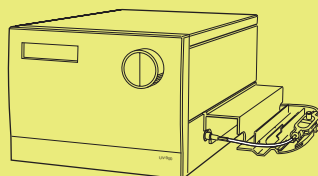

- 1 **Switch on the instrument** by means of the mains switch on the back panel.

Calibrating
Please wait...

- 2 **Switch on the UV-lamp** by selecting **Run**.

End
Press OK to **Run**

- 3 **Select wavelength** for one, two or three wavelengths.

Set Wavelength
(215, 254, 280nm)

Set Wavelength I1
(215nm)

Set Wavelength I2
(254nm)

Set Wavelength I3
(280nm)

- 4 **Autozero** the instrument by pressing the OK-button.

Autozero

- 5 The **main operating menu** (RUN-menu) is shown.

I1(215)	1.123
I2(254)	0.02345

