
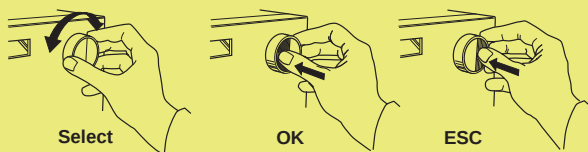


Short instructions Pump P-900

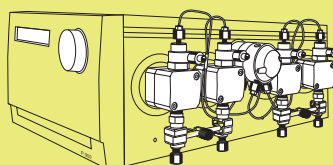

1 **Switch on the instrument** by means of the mains switch on the rear panel.

Selftest
Please wait...

2 Press **OK** to start synchronisation.

Pump synchronization
Press OK to start

3 The **main operating menu** is shown.

End 0.00ml/min
0.00MPa 0.0%B

4 Set the flow rate.

Set Flow Rate
(0.00ml/min) 2.5

5 Start the pump by setting it in **RUN**-mode.

End 2.50ml/min
Run

6 The **main operating menu** is shown.

Run 2.50ml/min
2.0MPa 0.0%B

7 Start a gradient by setting the **Gradient Length** and **Gradient Target**.

Set Gradient Length
(0.00min) 8.0

Set Gradient Target
(0.0%B) 10.0

8 Stop the pump by setting it in **END**-mode.

Run 2.50ml/min
End Hold Pause

