
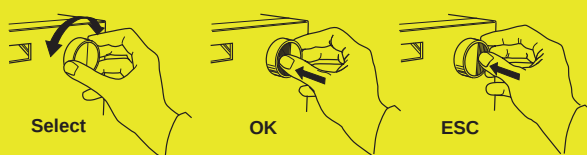


Short instructions Pump P-920

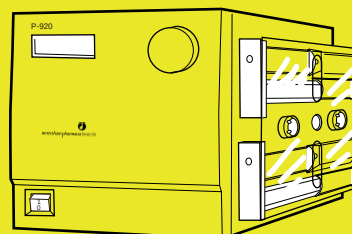

- 1 Switch on the module with the mains switch to the left on the front panel.
- 2 A self-test is performed. The version number is shown for two seconds.
- 3 The main operating menu is shown.
- 4 Set the flow rate.
- 5 Start the pump by setting it in **RUN** mode.
- 6 The main operating menu is shown.
- 7 Set concentration and gradient.
- 8 Start a gradient by setting the **Gradient Length** and **Gradient Target**.
- 9 Stop the pump by setting it in **END** mode.

Selftest
Please wait...

Pump P-920
<version no.>

End 0.00ml/min
0.00MPa 0.0%B

Set Flow Rate
(0.00ml/min) 2.5

End 2.50ml/min
Run

Run 2.50ml/min
2.0MPa 0.0%B

Set Conc./Gradient
(0.0%B) 10.0

Set Gradient Length
(0.00min) 8.0

Set Gradient Target
(0.0%B) 10.0

End 2.50ml/min
Run

Menu overview Pump P-920

