


Mounting kit for Pump P-950/P-960


Instruction

CAUTION! Before installing Pump P-950 or Pump P-960 in an ÄKTA™ design system, contact your local Amersham Biosciences service organisation to make sure that the system is equipped with the correct electrical components.


Installation

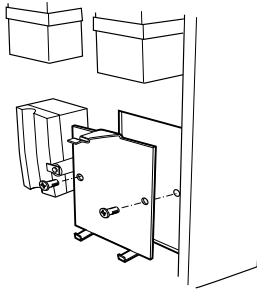


1. Attach the mounting plate to the lower right-hand corner at the inside of the system door, using the two flat-headed screws.

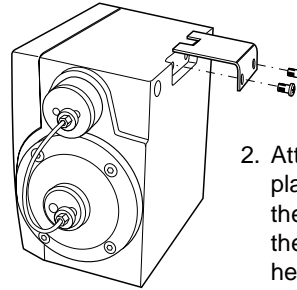

2. Attach the plate bracket to the pump using the two fillister-head screws.

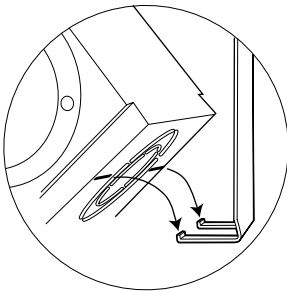


3. Fit the two hooks on the mounting plate in the holes underneath the pump.

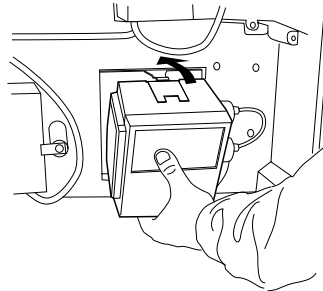

4. Push the pump towards the mounting plate as the figure shows. A click sound will indicate that the pump is locked in position.

Removing the pump

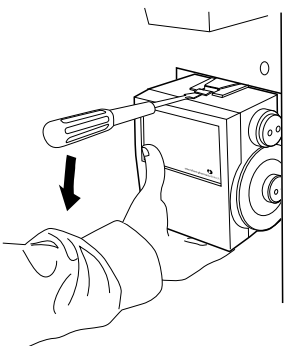

1. Release the locking plate with a screw driver while holding the pump.
2. Remove the pump.



WARNING! PINCH HAZARD
Do not grip the pump with the fingers at the rear of the pump.

Terms and Conditions of Sale

Unless otherwise agreed in writing, all goods and services are sold subject to the terms and conditions of sale of the company within the Amersham Biosciences group which supplies them. A copy of these terms and conditions is available on request.

Should you have any comments on this product, we will be pleased to receive them at:

Amersham Biosciences AB

SE-751 84 Uppsala
Sweden

Trademarks

ÅKTA is a trademark of Amersham Biosciences Limited.

Amersham and Amersham Biosciences are trademarks of Amersham plc.

© Copyright Amersham Biosciences AB 2003

- All rights reserved

Office Addresses

Amersham Biosciences AB

SE-751 84 Uppsala
Sweden

Amersham Biosciences UK Limited

Amersham Place
Little Chalfont
Buckinghamshire
England HP7 9NA

Amersham Biosciences Corp.

800 Centennial Avenue
PO Box 1327
Piscataway NJ 08855
USA

Amersham Biosciences Europe GmbH

Munzinger Strasse 9
D-79111 Freiburg
Germany

Amersham Biosciences K.K.

Sanken Building
3-25-1 Hyakunincho, Shinjuku-ku
Tokyo 169-0073
Japan

Amersham Biosciences Limited

13/E, Tower I
Ever Gain Plaza
88 Container Port Road
Kwai Chung, New Territories
Hong Kong

www.amershambiosciences.com