


# UNICORN 3.2 vs 3.1

---

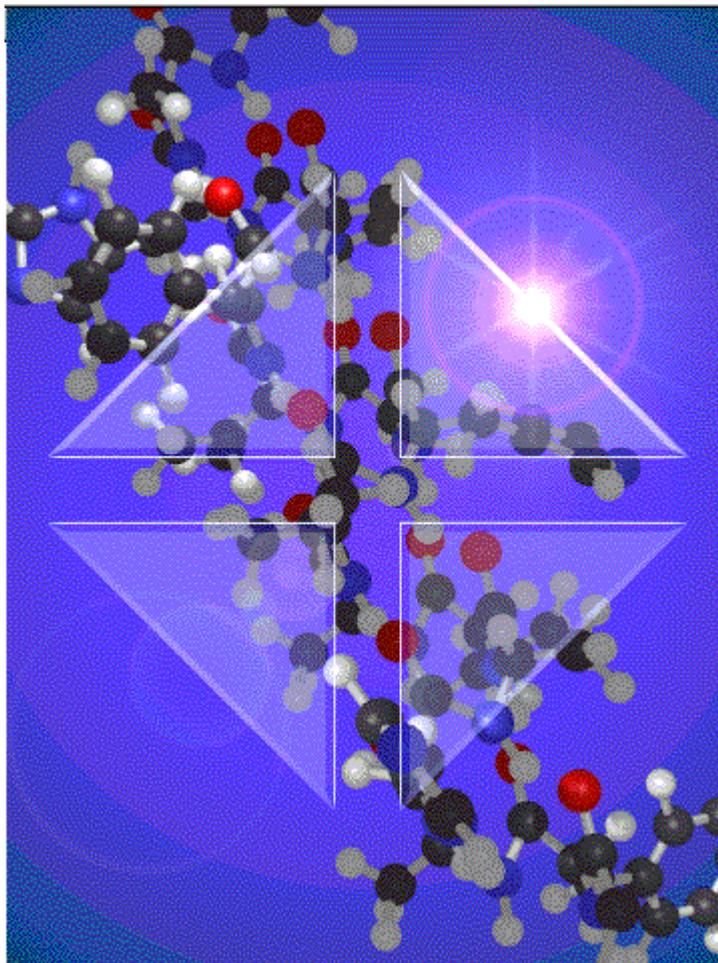

## Introduction

This document describes the implemented changes and improvements in UNICORN™ 3.2 compared to version 3.1.

**Note:** In some cases an updated strategy is needed in order for the new function to be available.

## Fraction collector

- UNICORN 3.2 supports the Frac-950 fraction collector.

## Software improvements

- The unicorn.ini file has been adjusted to avoid problems with the automatic evaluation of the installation test.
- The system will now update the accumulated volume for runs exceeding 20000 ml.
- FPLC™ strategies can now contain both **Watch\_Off** and **Hold\_Until** instructions.
- Gradients longer than 9999 ml are no longer accepted to avoid problems with pump capacity.
- Problems with the Evaluation module terminating when Oligo results are opened have been corrected.
- The **Customised Report** formats have been updated so the procedure **Quant\_Sample** can be used without causing the Evaluation module to terminate.
- The Edit system dialog has been adjusted so that the correct strategy is selected when the strategy has a name that is a sub name of another strategy.
- The Method Notes have been expanded to accept up to 20000 characters.

## Important Information

ÄKTA, UNICORN, FPLC and Dropdesign are trademarks of Amersham Biosciences Limited.

Amersham and Amersham Biosciences are trademarks of Amersham plc.

Windows NT and Windows 2000 are trademarks of Microsoft Corporation.

Compaq is a trademark of Compaq Computer Corporation. Novell and NetWare are registered trademarks of Novell Inc.

© Amersham Biosciences AB 2002 – All rights reserved

All goods and services are sold subject to the terms and conditions of sale of the company within the Amersham Biosciences group which supplies them. A copy of these terms and conditions is available on request.

### **Amersham Biosciences AB**

Björkgatan 30  
SE-751 84 Uppsala  
Sweden

### **Amersham Biosciences Europe GmbH**

Munzinger Strasse 9  
D-79111 Freiburg  
Germany

### **Amersham Biosciences China Limited**

13/F., Tower I  
Ever Gain Plaza  
88 Container Port Road  
Kwai Chung, New Territories  
Hong Kong

### **Amersham Biosciences UK Limited**

Amersham Place  
Little Chalfont  
Buckinghamshire  
England HP7 9NA

### **Amersham Biosciences Corp**

800 Centennial Avenue  
P.O. Box 1327  
Piscataway, NJ 08855  
USA

### **Amersham Biosciences K.K.**

Sanken Building, 3-25-1  
Hyakunincho, Shinjuku-ku  
Tokyo 169-0073  
Japan.

**[www.amershambiosciences.com](http://www.amershambiosciences.com)**

